

Collection Bucket Decontamination Protocol:

Materials:

- Collection buckets
 - Bleach
 - Tap water
 - Deionized [DI] water
 - Sponge
-

Protocol

1. Scrub your first bucket [B1] and lid w/ 10% bleach solution in crevices using a sponge
2. Fill B1 with 10% bleach, ½ way, seal with lid
3. Shake B1 for 15 seconds, then let sit for 5 minutes
4. Scrub bucket 2 [B2] while waiting for B1,
5. Pour bleach solution from B1 to B2, then shake and let sit
6. Rinse B1 with tap water
7. Fill B1 ½ way with tap water and shake (2x)
8. Rinse B1 with DI water
9. Fill B1 ½ way with DI water and shake (2x)
10. Repeat 4-10 until the desired number of buckets are clean